## **Gebrauchsanweisung Instruction manual Manual de instrucciones**

### **Taumel-Rollenmischer mit 5 Rollen Swivel roller mixers with 5 rollers Agitadores orbitales con 5 rollos „Swap“-System**

#### **RM5W-80V**

- herausnehmbare Rollen zur Schaffung von Liegefläche für Gefäße mit einem Volumen von bis zu 5 Liter
- Umbau ohne Werkzeug
- Drehzahl regelbar von 10 bis 80 U/min.
  
- removable rollers for vessels with larger diameter with a volume of up to 5 litres
- tool-free change of rollers
- speed infinitely adjustable from 10 to 80 rpm
  
- rollos extraíbles para vasos con mayor diámetro con un volumen de hasta 5 litros
- cambio de rollos sin herramientas
- velocidad ajustable continuamente de 10 a 80 r.p.m.

## Inhalt / Content

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## 2. General information

The roller mixers are designed in accordance with Safety Class 1. They are produced and tested in accordance with DIN EN 61010. According to these regulations, the units are designed to meet the requirements for safe and correct operations. To ensure the proper safety and operational functions of the instrument, the user should follow the instructions and safety guidelines in this manual.

### 2.1. Unpacking the instrument

Unpack the instrument carefully and check to see that it is not damaged. It is important that any damage incurred during transport, such as indented sides of the carton, chemical impairment of the carton, water stains or other physical impacts which might have damaged the content, be recognized at the time of unpacking. Notify your carrier or forwarding agent immediately in case of such damage. You may unpack the carton carefully if there is no visible damage of the packing material.

#### Note:

If the instrument is not damaged and all parts are complete, you may start to operate the device after reading the instruction manual.

#### Important note:

Read this instruction manual carefully before operating the instrument. Should there be any further questions after reading the manual concerning the setup, operation or warranty, please contact:

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

**Important note:**


Please store the manual in a place easily accessible for every user.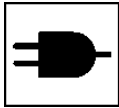

A EURO AC plug (DIN 49441 CEE 7/VII10/ 16 A 250 V) is normally shipped with our instruments. North America instruments are shipped with standard US plugs (NEMA Pub. No. WDI1961 ASA C 73.1. 1961 page 8 15A 125V).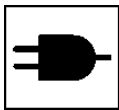

When operating the instruments in countries with different AC plug systems, use an approved adapter or have a qualified electrician replace the AC plug with an approved model suitable for the country of operation. Ensure that your local supply voltage matches the indication on the instrument's specification plate!

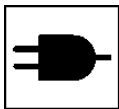The instrument is earthed. When replacing the original AC plug, ensure that the earth conductor is connected to the new plug!

### 3. Safety instructions

#### 3.1 Warning icons on the instrument

**Warning against general danger:**

This symbol indicates that it is imperative to read and understand the instruction manual prior to operating the instrument. Please highlight points which require special attention in your field of application so they are not overlooked. Disregard of warnings may result in impairment of serviceability as well as in physical harm to the user.

### 3.2 Warning icons in the instruction manual

This symbol distinguishes notices which have to be strictly observed by the user. Disregard of warnings may result in the impairment of serviceability as well as in physical harm to the user.

This symbol distinguishes notices which have to be strictly observed by the user to ensure safe operation of the unit.

Note regarding repair or maintenance

Note regarding mains voltage

### 3.3 General safety instructions

**Caution:**  
Please comply with all safety and accident-prevention regulations applicable to laboratory work.

**Caution:**  
Please do not use this instrument in insecure environment especially not in explosive ambience. Danger of life!

**Caution:**  
Instructed users only may operate the instrument.

**Caution:**  
When connecting the instrument to an AC power outlet, ensure that your local supply voltage corresponds to the specifications indicated on the instrument.


**Caution:**  
Please act with caution when working in the vicinity of flammable or explosive substances and observe safety data sheets. This instrument is

not explosion-proof.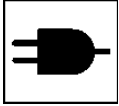

The On/Off switch does not disconnect the instrument from the power source. Remove the plug from the AC power outlet to disconnect the instrument from the mains supply entirely.


**Note:**

Make sure that all safety information on the instrument is clearly visible during operation.

**Attention:**

Do not open the instrument. Repairs are only to be carried out by trained service technicians.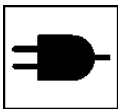

Connect the instrument to an earthed AC power outlet only.

Make sure the unit is standing on a solid and horizontal surface.

#### 4. Intended use

The roller mixers are mainly used to mix hydrous solutions not containing inflammable parts in glass or plastic vessels. They are designed for the use in chemical and biological laboratories of industrial enterprises, universities and pharmacies.

To ensure maximum service life, observe the specified ambient conditions (temperature and humidity) and ensure that the instrument is not exposed to a corrosive atmosphere.

## 5. Setup

**Caution:**

Please do not use this instrument in insecure environment especially not in explosive ambience. Danger of life!

**Caution:**

When connecting the instrument to an AC power outlet, ensure that your local supply voltage corresponds to the specifications indicated on the instrument.

Rocker switch (1)

Speed control button (2)

### 5.1 Operation "Mixing"

Five synchronised rotating PVC rollers form a homogeneous reel roller bearing for the use of closed test tubes. The tubes are both, simultaneously rotated along their longitudinal axis and rocked back and forth on their bearings.

Start the mixing function by means of the rocker switch (1) ON/OFF. Set speed by means of the speed control button (2).

### 5.2 Cleaning and maintenance

The housing and the 5 PVC rollers are acid-resistance covered for long lasting use and easy cleaning with warm water and dish washing liquid. The instrument is maintenance-free.

It must be opened and repaired only by authorized service personnel. Any work on the electronics in the unit should only be carried out by knowledgeable, trained personnel.

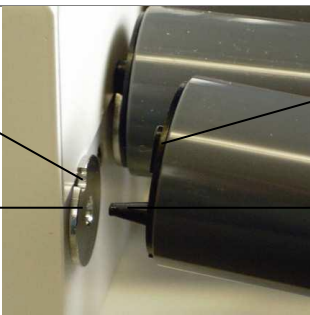
**Caution:**


Unplug the AC power cord before opening the housing.

### 5.3 Removing the rollers

 <p>Knurled head screw</p>	<p>a) Switch off instrument</p> <p>b) Loosen knurled head screw by 4 rotations to the left.</p> <p><b>Do not</b> remove the screw completely!</p>
	<p>c) The coverage opens</p>
	<p>d) The exchange of the rolls may be carried out optionally. Lift the rollers carefully out of the bearing bar. The roller now can be removed axially.</p>
	<p>e) When replacing the rollers please follow steps a) to c)</p> <p>f) Then push the conical part of the roller (1) into the borehole of the drive disc (2). Then rotate the roller axial until the drive piece (3) clicks into the slot of the drive disc (4)</p>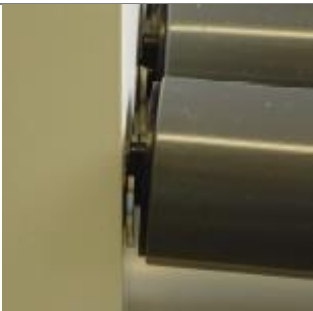

			<p>g) Once the rollers are replaced...</p>
			<p>h) .... close the coverage with the knurled head screw. Take care that the coverage is completely closed.</p>

## 6. Dismantling, transport and storage

### 6.1 Dismantling

1. Switch off the instrument by means of the rocker switch.
2. Disconnect the instrument from the mains.
3. Now you can remove the instrument from the working area.

### 6.2 Transport and Storage

Place the instrument and its parts in its original packaging or in a suitable container to protect it during transport. Close the packaging with adhesive tape.

In case you do not use the original packaging, please mark the box with the following notes:

- Glass symbol (Fragile! Handle with care!)
- Umbrella (Keep dry!)
- Content (List of content)

Please observe the following conditions of surrounding:

max. ambient temperature	+5-40 °C
max. ambient humidity	80 %

## 7. Disposal

Damaged and/or waste electric or electronic equipment have to be disposed of at the designated Recycling Depot.

## 8. Warranty and Liability

Paul Marienfeld GmbH & Co. KG provides a guarantee for a period of 24 months from date of delivery for the instrument purchased by you. This warranty applies only to defects in material or workmanship. There is no warranty for defects or malfunctions caused by neglect, improper use or improper service.

In case of defects in material or workmanship, the instrument will be repaired or replaced. In case of replacement, the warranty for the new unit will persist for the remaining period of the primary warranty period. Charges for returning the instrument have to be borne by the customer.

Any claims regarding guarantee and liability are limited to professional and proper use of the instrument and observance of the general safety instructions and this instruction manual.

Paul Marienfeld GmbH & Co. KG is not liable for any material or personal damages caused by improper use and does not accept liability for consequential damages. The warranty expires in case of removal or alteration of the identification label and/or the serial number of the instrument.

## 9. Technical Data

Mains	115 resp. 230 Volt, 50-60 Hz (see rear panel)
Power	25 Watt
Speed	10-80 U/min.
Length of rollers	330 mm
Diameter of rollers	32 mm
Ø distance from roller to roller	36.5 mm
Dimensions	436x177x105 mm
Weight	3.8 kg
Max. load	1.5 kg
Max. ambient temperature	+5 °C bis +40 °C
Max. operating period	100 %

## 10. Declaration of conformity

We hereby declare that the following products

Swivel roller mixer RM5W-80V, 230 V (Cat. no. 72 040 28) or  
Swivel roller mixer RM5W-80V, 115 V (Cat. no. 72 140 28)

correspond to the regulations

2006/95/EG  
2004/108/EG  
2011/65/EU und  
2006/42/EG

and are in accordance with the standards or standardized documents:

DIN EN 61010-1  
DIN EN 61010-2-051  
DIN EN 61326-1

In the case of a modification of the unit which has not been agreed on with us, this declaration becomes null and void and the warranty expires.

Paul Marienfeld GmbH & Co. KG  
Gerd Marienfeld, President<sup>1</sup>

<sup>1</sup> The above declaration has been generated automatically and is valid without signature

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