## **Gebrauchsanweisung Instruction manual Manual de instrucciones**

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## 2. General information

The roller mixers are designed in accordance with Safety Class III (EN 6114). They are produced and tested in accordance with DIN EN 61010. According to these regulations, the units are designed to meet the requirements for safe and correct operations. To ensure the proper safety and operational functions of the instrument, the user should follow the instructions and safety guidelines in this manual.

### 2.1. Unpacking the instrument

Unpack the instrument carefully and check to see that it is not damaged. It is important that any damage incurred during transport, such as indented sides of the carton, chemical impairment of the carton, water stains or other physical impacts which might have damaged the content, be recognized at the time of unpacking. Notify your carrier or forwarding agent immediately in case of such damage. You may unpack the carton carefully if there is no visible damage of the packing material.

#### Note:

If the instrument is not damaged and all parts are complete, you may start to operate the device after reading the instruction manual.

#### Important note:

Read this instruction manual carefully and make sure everything is understood before operating the instrument. Should there be any further questions after reading the manual concerning the setup, operation or warranty, please contact:

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
**Important note:**

Please store the manual in a place easily accessible for every user.

**3. Safety instructions****3.1 Warning icons on the instrument****Warning against general danger:**

This symbol indicates that it is imperative to read and understand the instruction manual prior to operating the instrument. Please highlight points which require special attention in your field of application so they are not overlooked. Disregard of warnings may result in impairment of serviceability as well as in physical harm to the user.

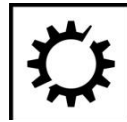**3.2 Warning icons in the instruction manual****3.2.1 Explanation of the symbols****Hazard point****Bio hazard****Risk of fire****Explosion****Mechanical hazard****Material damage****3.2.2 Explanation of the signal words****Danger**

Signifies a dangerous situation. Non-observance can lead to severe injuries or death.

**Warning**

Indicates a possibly dangerous situation. Non-observance can lead to severe injuries or death.

**Caution**

Indicates a possibly dangerous situation. Non-observance may lead to light or moderate injuries.

**Notice** Please observe the notice. Non-observance may lead to material damage.

### 3.3 General safety warnings and instructions

#### **Danger: Risk of explosion.**

- Please do not use this instrument in insecure environment, especially not in explosive ambience. Danger to life!
- Do not use this instrument next to highly flammable or explosive substances. This instrument is not explosion-proof!
- Do not use this device for processing any substances which could generate an explosive atmosphere.
- Do not use this device to process any explosive or highly reactive substances.

#### **Warning: Damages to health due to infectious liquids and pathogenic germs**

- When handling infectious liquids and pathogenic germs, observe the national regulations, the biological security level of your laboratory, the material safety data sheets and the manufacturer's application notes.
- Wear personal protective equipment.
- For comprehensive regulations about handling germs and biological material of the risk group II or higher, please refer to the "Laboratory Biosafety Manual" in its respectively current valid version from the World Health Organisation.

#### **Warning: Risk of fire**

Act with caution handling highly flammable liquids. Please observe the safety data sheets. This instrument is not explosion-proof!

#### **Warning: Damage to health due to contaminated device and accessories**

In the following cases sample material can be released:

- improperly sealed tubes
- instable tubes
- high vapor pressure of the content so that the seal of the tubes can spring open
- damaged sealing
- smashed glass tubes
- Only mix in closed tubes.
- Observe the nationally prescribed safety environment when working with hazardous, toxic and pathogenic samples. Pay particular attention to personal protective equipment (gloves, clothing, goggles etc.), extraction and the safety class of the lab.

- Decontaminate the device and the accessories before storage and shipping.

**Caution: Poor safety due to incorrect accessories and spare parts.**

The use of accessories and spare parts other than recommended by Paul Marienfeld GmbH & Co. KG may impair the safety, function and precision of the device. Paul Marienfeld GmbH & Co. KG cannot be held liable or accept any liability for damage resulting from the use of incorrect or non-recommended accessories and spare parts, or from the improper use of such equipment.

- Only use accessories and spare parts recommended by Paul Marienfeld GmbH & Co. KG.

**Caution: Crush hazard due to moving parts.**

- Do not replace any consumables as long as the device is running.
- Do not open the coverage whilst the device is running.
- Only switch-on the device with properly mounted rolls and completely closed coverage.
- Do not place hands or objects into the area under the rollers while the machine is running.

**Notice: Damage to electronic components due to condensation.**

Condensate can form in the device after it has been moved from a cool environment to a warmer one.

- Wait at least for 3 hours after installing the device before connecting it to mains.

**Notice: Damage to electronic components due to incorrect supply voltage.**

Incorrect supply voltage may destroy electrical components.

- Only use a power supply with a 24 V DC, min. 600 mA output

**Notice: Damage from the use of aggressive chemicals.**

- Do not use any aggressive chemicals on the device or its accessories such as strong or weak bases, acids, acetone formaldehyde, halogenated hydrocarbons or phenol.
- If the device has been contaminated by aggressive chemicals, immediately clean it by means of a mild cleaning agent.

#### 4. Intended use

The roller mixers are mainly used to mix hydrous solutions not containing inflammable parts in glass or plastic vessels. They are designed for the use in chemical and biological laboratories of industrial enterprises, universities and pharmacies.

To ensure maximum service life, observe the specified ambient conditions (temperature and humidity) and ensure that the instrument is not exposed to a corrosive atmosphere.

#### 5. Setup

##### Caution:

Please do not use this instrument in insecure environment especially not in explosive ambience. Danger of life!

##### Notice:

The roller mixer is powered by a 24 Volt DC power supply with an angled hollow plug 5.5/2.1 mm.  Make sure to use a suitable power supply only.

The input socket is located on the bottom of the housing, the On/Off switch (1) on the side of the housing. The speed is set with the speed control button (2).

Rocker switch  
(1)

Speed control  
button (2)

### 5.1 Operation „Mixing“

Five resp. ten synchronised rotating PVC rollers form a homogeneous reel roller bearing for the use of closed test tubes. The tubes are both, simultaneously, rotated along their longitudinal axis and rocked back and forth on their bearings.

Start the mixing function by means of the rocker switch (1) On/Off. Set speed between 2 U/min. and 70 U/min. (steplessly) by means of the speed control button (2).

### 5.2 Cleaning and maintenance


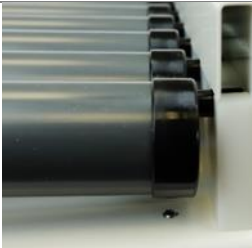
The housing and the 5 resp. 10 PVC rollers are acid-resistance covered for long lasting use and easy cleaning with warm water and dish washing liquid.

The instrument is maintenance-free. It must be opened and repaired only by authorized service personnel. Any work on the electronics in the unit should only be carried out by knowledgeable, trained personnel.

### 5.3 Removing and inserting of individual rolls

1		<p>a. Switch off the device. b. Loosen knurled head screws by 4 rotations to the left.</p>
2		<p>c. The coverage opens.</p>
3		<p>d. The exchange of rolls may be carried out optionally. Lift the rolls carefully out of the bearing bar. The roll now can be removed axially. Thereafter close the coverage again.</p>

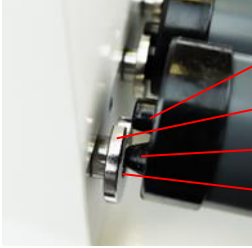4		<p>e. When replacing the rolls please follow steps a to c.</p> <p>f. Then push the conical part of the roll (1) into the borehole of the drive disk (2). Now rotate the roll axial until the drive piece (3) clicks into the slot of the drive disk (4).</p>
5		<p>g. Take care that the rolls are completely inserted into the drive disk.</p>
6		<p>h. Close the coverage with the knurled head screw. Take care that the coverage is completely closed.</p>

## 6. Dismantling, transport and storage

### 6.1 Dismantling

1. Switch off the instrument by means of the rocker switch.
2. Disconnect the instrument from the mains.
3. Now you can remove the instrument from the working area.

### 6.2 Transport and storage

Place the instrument and its parts in its original packaging or in a suitable container. Protect each module by means of bubble wrap against mechanical shocks from the outside. When packing the individually packed modules in one carton, make sure there are sufficient gaps between them and fill out the gaps with insulating material. In case you do not use the original packaging, please mark the box with the following notes:

- Glass symbol (Fragile! Handle with care!)
- Umbrella (Keep dry!)
- Content (List of content)

For proper storage please take note of the following instruction:

1. For storage of the device, separate the individual modules.
2. Pack each module into a separate, air-tight plastic bag.
3. Please observe the following conditions of surrounding:
 

max. ambient temperature	+5-40 °C
max. ambient humidity	80 %

## 7. Disposal

Damaged and/or disposed electric or electronic equipment has to be disposed of at the designated Recycling Depot. Also the packing material should be disposed of environment-friendly (material separation).

## 8. Warranty and Liability

Paul Marienfeld GmbH & Co. KG provides a guarantee for a period of 24 months from date of delivery for the instrument purchased by you. This warranty applies only to defects in material or workmanship. There is no warranty for defects or malfunctions caused by neglect, improper use or improper service.

In case of defects in material or workmanship, the instrument will be repaired or replaced. In case of replacement, the warranty for the new unit will persist for the remaining period of primary warranty period. Charges for returning the instrument must be borne by the customer.

Any claims regarding guarantee and liability are limited to professional and proper use of the instrument and observance of the general safety instructions and this instruction manual.

Paul Marienfeld GmbH & Co. KG is not liable for any material or personal damages caused by improper use and does not accept liability for consequential damages. The warranty expires in case of removal or alteration of the identification label and/or the serial number of the instrument.

## 9. Technical Data

	<b>RM5</b>	<b>RM10</b>
Mains	24 V DC	24 V DC
Power	15 Watt	15 Watt
Speed	10-70 U/min.	10-70 U/min.
Length of rollers	330 mm	330 mm
Diameter of rollers	32 mm	32 mm
Distance from roller to roller	36 mm	36 mm
Dimensions	480x180x113 cm	480x360x113 cm
Weight	3.7 kg	6.5 kg
Max. load	5.0 kg	5.0 kg
Max. ambient temperature	+5 °C bis +40 °C	+5 °C bis +40 °C
Max. operating temperature	100 %	100 %
Max. ambient relative humidity	80 %	80 %

## 10. EC-Declaration of conformity in accordance with EEC machine directive 2006/42/EG appendix II 1A

We declare under our sole responsibility that the following product

Roller mixer RM5/RM10 | 24VDC | 15 W

corresponds to the regulations and conforms with the standards or standardized documents.

The agreement with further valid guidelines/regulations for the product is explained:

EMV – Directive 2014/30/EG

RoHS – Directive 2011/65/EU

Applied harmonized standards in particular:

DIN EN 61326-1: 2013-07

DIN EN 61010-1: 2011-07

DIN EN ISO 12100: 2011-03

DIN EN 61010-2-051: 2016-02

In case of a modification of the unit which has not been agreed on with us, this declaration becomes null and void and the warranty expires.

Paul Marienfeld GmbH & Co. KG

Harry Marienfeld, Director <sup>1</sup>

<sup>1</sup> The above declaration has been generated automatically and is valid without signature

## User Instructions Power supply

### Proper use

This AC/DC adapter is designed as external power supply for electrical equipment. It may only be used indoors.

### Installation

Please read and follow the instructions before first use.

Before first use please check the mains voltage to match the input voltage specified on the power supply. Then check the output voltage, its polarity, the output current and power to match the requirements of the equipment to be connected with.

The connection of the plug must be stable; twisting or bending of the cable should be avoided. Provide adequate ventilation during operation. The adapter should not be covered or used near the heating sources and may not be exposed to direct sunlight.

If your device is not used for a long time we recommend to disconnect the adapter from main voltage in order to avoid any damage due to overvoltage for example (spikes, lightning).

### Maintenance

For cleaning you may not use cleaning agents. Wipe with dry cloth. Disconnect the adapter from the mains before cleaning.

### Disposal

Electronic devices are hazardous waste and do not belong to the normal house waste. If the device reached end of its life dispose it according to the applicable legal provisions of the municipal collection points. Disposing in the domestic waste is not permitted.

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MEAN WELL ENTERPRISE Co., Ltd.
Schukat electronic Vertriebs GmbH