Trockenbank für Objektträger
Drying Bench for Microscope Slides
Plataforma mecánica para secar portaobjetos

H 15/30

Inhalt

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Content

1. General information	13
1.1 Unpacking the instrument	13
1.2 Scope of delivery	13
2. Safety instructions	14-15
2.1 Description of safety symbols in the instruction manual	14
2.2 Description of safety symbols on the instrument	14
2.3 General safety instructions	15
3. Intended use	16
4. Setting up the instrument	16
5. Operation	16, 17
6. Maintenance and service	17-18
6.1 Service and repair	17, 18
6.2 Cleaning	18
7. Disassembly, packing and storage	18-19
7.1 Disassembly	18
7.2 Packing and storage	18, 19
8. Repair	19
9. Technical data	19
10. Warranty and liability	19
11. Liability exclusion	19
12. Disposal	19

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1. General information

1.1 Unpacking the instrument

Our slide drying bench H 15/30 has been manufactured and tested according to DIN EN 61010. According to these regulations, the units are designed to meet the requirements for safe and correct operations. To ensure the proper safety and operational functions of the instrument, the user should follow the instructions and safety guidelines in this manual.

The instrument comes with a SCHUKO plug (DIN 49441, 16 A, 250 V). When operating the instrument in countries with different AC plug systems, use an approved adapter or have a qualified electrician replace the AC plug with an approved model suitable for the country of operation.

Unpack the instrument carefully and check to see that it is not damaged. It is important that any damage incurred during transport be recognized at the time of unpacking. Notify your carrier or forwarding agent immediately in case of such damage.

1.2 Scope of delivery

Your delivery should contain the following:

- 1 Slide drying bench H 15/30
- 1 Slide support
- 1 Instruction manual

Note:

If the instrument is not damaged and all parts are complete, you may start to operate the device after reading the instruction manual.

Important note:

Read this instruction manual carefully before operating the instrument. Please store the manual in a place easily accessible for every user. Should there be any further questions after reading the manual concerning the setup, operation or warranty, please contact:

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2. Safety instructions

2.1 Description of safety symbols in the instruction manual

This symbol distinguishes notices which have to be strictly observed by the user. Disregard of warnings may result in the impairment of serviceability as well as in physical harm to the user.

This symbol indicates proscriptions, which have to be strictly observed by the user. Disregard of the proscriptions may result in the impairment of serviceability as well as in physical harm to the user.

This symbol distinguishes notices which have to be strictly observed by the user to ensure safe operation of the unit.

Caution
Danger for fire or explosion!

Note regarding repair or maintenance

Note regarding mains cable connection

2.2 Description of safety symbols on the instrument

Warning against general danger: This symbol indicates that it is imperative to read and understand the instruction manual prior to operating the instrument. Please highlight points which require special attention in your field of application so they are not overlooked. Disregard of warnings may result in impairment of serviceability as well as in physical harm to the user.

Warning against hot surfaces: This symbol indicates hot surfaces on the device. It refers to the hotplate of the instrument which can reach up to 115 °C during operation. There is danger of burns for the user and danger of ignition for flammable substances which come into contact with the heating surface or which are in the immediate proximity.

2.3 General safety instructions

Caution:

Please comply with all safety and accident-prevention regulations applicable to laboratory work.

Caution:

Please do not use this instrument in insecure environment especially not in explosive ambience. Danger of life!

Caution:

Instructed users only may operate the instrument.

Caution:

When connecting the instrument to an AC power outlet, ensure that your local supply voltage corresponds to the specifications indicated on the instrument.

Caution:

Please act with caution when working in the vicinity of flammable or explosive substances and observe safety data sheets. This instrument is not explosion-proof.

The On/Off switch does not disconnect the instrument from the power source. Remove the plug from the AC power outlet to disconnect the instrument from the mains supply entirely.

Note:

Make sure that all safety information on the instrument is clearly visible during operation.

Attention:

Do not open the instrument. Repairs are only to be carried out by trained service technicians.

Make sure the unit is standing on a solid and horizontal surface.

3. Intended use

The slide drying bench H 15/30 is designed to dry up to 48 microscope slides in various positions which are described in chapter 5.

4. Setting up the instrument

The instrument is designed to be safe under the following conditions

- indoor use
- temperatures between 5 °C and 40 °C
- max. relative humidity 80 % with no condensation inside and outside of the housing
- mains supply voltage fluctuations up to ± 10 % of the nominal voltage

The instrument has to be used with a power cable with protective conductor and rubber connector.

When connecting the instrument to an AC power outlet, make sure that your local supply voltage matches the indication on the instrument.

Ensure only the correct rated mains input fuses are fitted – see chapter 9 “Technical data”

Install the instrument on a clean, dry- non-combustible, solid work surface with a tleast 300 mm distance from other equipment.

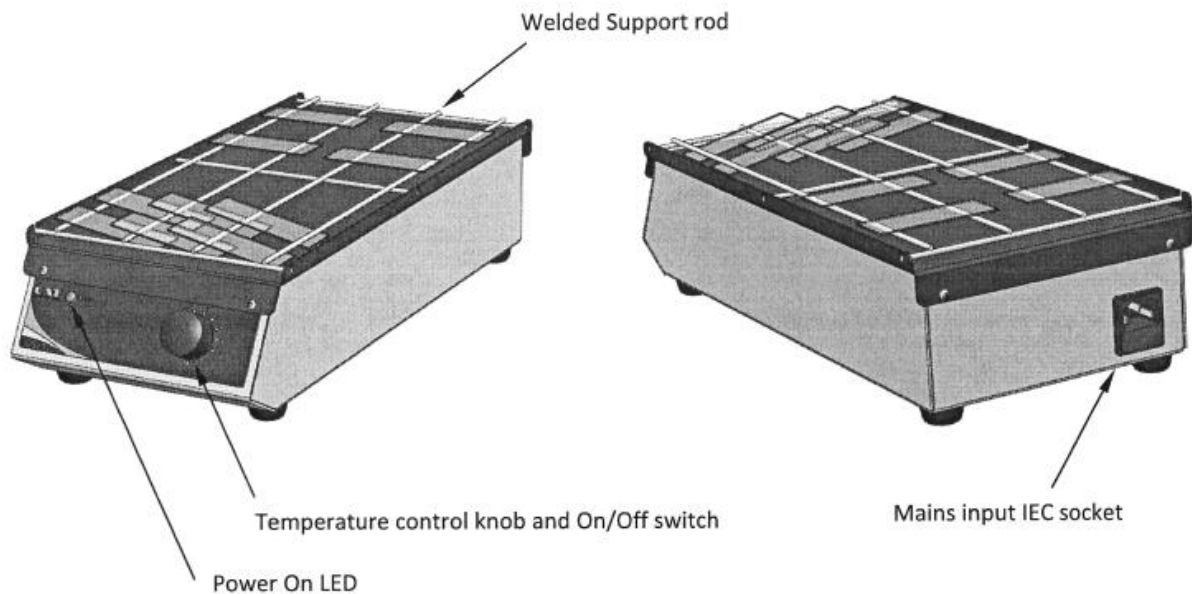
Important:

Failure to use this equipment in accordance with the manufacturer's operation manual may compromise your basic safety protection afforded by this equipment and may invalidate the terms of the warranty/guarantee.


5. Operation

Place the slides to be dried on the bench surface. The slides can be arranged on the bench in a number of ways as illustrated following herewith, with or without the welded support.

- Connect the device to the mains power ensuring the correct voltage is being applied.
- Set the temperature control knob to the desired temperature. The power on LED will light up.
- The max. temperature of the bench surface is 115 °C
- When slides are dry, switch off the unit by means of the regulator knob and disconnect from the mains electrical supply.
- Allow the slides to cool down before further processing

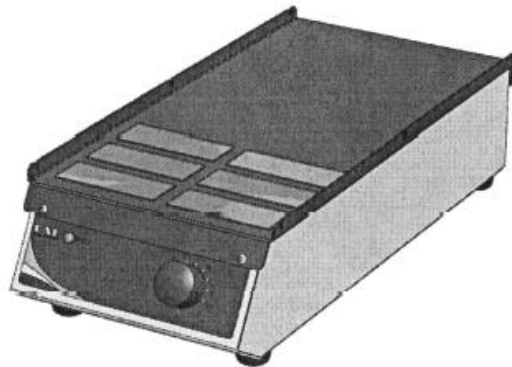Place the slides to be dried on the bench surface. The slides can be arranged on the bench in a number of ways as illustrated and without the welded support rod as well as shown in the following illustration.

6. Maintenance and service

6.1 Service and repair

Should you have any questions about installation, operation or maintenance, please contact the supplier of the device at the following address:

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**Caution:**

In case of malfunction do not attempt to repair the unit! There are no user-serviceable parts in this instrument. Any attempt by the user to repair the unit will immediately render the guarantee null and void.

**Caution:**

Do not open the device! Repair is only to be carried out by trained service technicians.

**Caution:**

In case of repair: If you are obliged to send the device back to the supplier, please make sure that the device has been adequately cleaned / decontaminated beforehand. The user is responsible for direct or indirect damage both corporal and material that may result from the media used.

6.2 Cleaning

The coated aluminium housing and polyamide plastic covers allow an easy cleaning of the instrument. The front panel is chemically stable and splash-proof. We advise to use a lint-free cloth and a mild soap solution (water and household detergent) for cleaning.

Please do not use steel-wool or aggressive detergent, otherwise damage to the lettering and surface cannot be avoided.

7. Disassembly, packing and storage

7.1 Disassembly

Switch off the device and disconnect it from the mains.

7.2 Packing and storage of the device

If possible, please use the original packing material for packing the instrument. Protect each module against shocks from outside with bubble wrap. If you pack individual modules in a box ensure adequate spacing between them and fill the spaces with insulation material. If you don't use the original packing, please mark the pack on the outside as follows:

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Observe the following notes for the correct storage of the instrument:

1. Separate the individual modules from each other for storing
2. Pack each module in a separate airtight plastic bag
3. Ensure the following ambient conditions:
 - max. ambient temperature: room temperature up to +40 °C
 - max. air humidity: 80 %

8. Repair

In case that the instrument came into contact with hazardous substances, please provide, when returning for repair, the following:

- precise information on the relevant medium
- protective measures to ensure the safety of our receiving and maintenance personnel
- the package marked as appropriate for hazardous materials

9. Technical data

Power supply	230 V AC or 115/100 V AC, 50/60 Hz
Max. power consumption	150 W
Fuse	230 V AC: IEC 60127-2 5x20 mm 2A T/250V 115 V AC: IEC 60127-2 5x20 mm 4A T/250 V
Heating element	silicone rubber mat heater
Ambient temperature range	5-40 °C
Max. humidity	80 % RH
Capacity	max. 48 microscope slides
Dimensions	395 x 186 x 105 mm
Weight	1.3 kg
Thermal insulation	Ceramic fibre sheet

10. Warranty and liability

Paul Marienfeld GmbH & Co. KG provides a guarantee for a period of 24 months from date of delivery for the instrument purchased by you. This warranty applies only to defects in material or workmanship. There is no warranty for defects or malfunctions caused by neglect, improper use or improper service.

In case of defects in material or workmanship, the instrument will be repaired or replaced. In case of replacement, the warranty for the new unit will persist for the remaining period of the primary warranty period. Charges for returning the instrument have to be borne by the customer.

Any claims regarding guarantee and liability are limited to professional and proper use of the instrument and observance of the general safety instructions and this instruction manual.

Paul Marienfeld GmbH & Co. KG is not liable for any material or personal damages caused by improper use and does not accept liability for consequential damages. The warranty expires in case of removal or alteration of the identification label and/or the serial number of the instrument.

11. Liability exclusion

Paul Marienfeld GmbH & Co. KG is not liable for any material or personal damages caused by improper use and does not accept liability for consequential damages. The warranty expires in case of removal or alteration of the identification label and/or the serial number of the instrument.

12. Disposal

Damaged and/or waste electric or electronic equipment has to be disposed of at the designated recycling depot.

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