Bedienungsanleitung
Instruction manual
Manual de instrucciones

Schüttelgerät S20
Shaker S20
Agitador S20

Inhalt / Content / Contenido

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Abmessungen: 130 x 132 x 40 mm
Gewicht: 0,140 kg


1. Scope of delivery

- 1 Shaker S20 (Drive unit without attachment)
Cat. no. 7204010 (230V/50-60Hz) or Cat. no. 7214010 (115V/50-60Hz)
- 1 Instruction manual


Attachments (have to be ordered separately):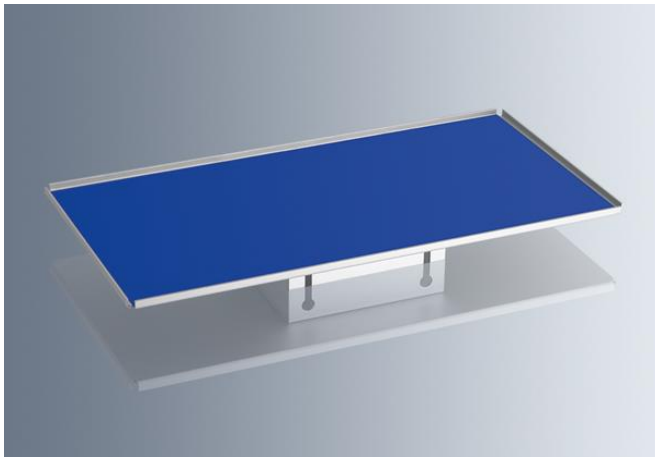

Shaking platform (Cat. No. 7204011):

This large platform (210x410 mm) has special adhesive mat which securely holds petri dishes, microtiter plates and other small beakers during shaking, without the need for fasteners.

Recommended max. shaking speed:
1100 rpm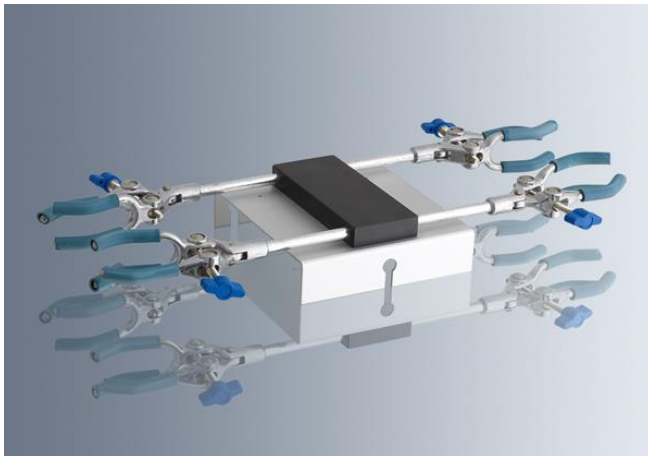

Four-flask shaker attachment (Cat. No. 7204012):

This attachment holds up to four Erlenmeyer-flasks (500 ml each) or round bottom flasks for slow to intensive, constant rotary motion. It is designed to provide a favourable centre of mass because of the lateral arrangements of the flasks.


Recommended max. shaking speed:
500 rpm

**Universal shaker attachment
(Cat. No. 7204013):**

This attachment comes with one fixed and one flexible elastic rubber roller to hold Erlenmeyer flasks, flat bottom flasks, round and square bottles and other various containers. All containers are held firmly in place for any constant, slow or fast speed setting.


Recommended max. shaking speed:
500 rpm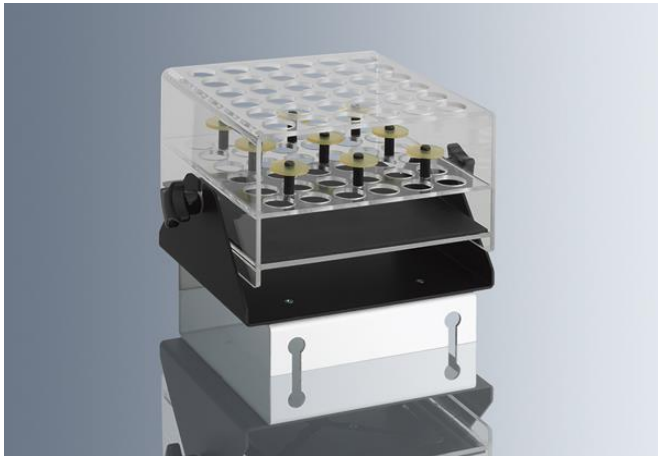

**Test tube attachment
(Cat. No. 7204014):**

This attachment provides for extremely slow to intensive, constant rotary motion for shaking up to 36 test tubes of up to 16 mm diameter. The two screws at the side allow for setting the angle of the test tubes from vertical to approximately 45° inclination.

Recommended max. shaking speed:
800 rpm

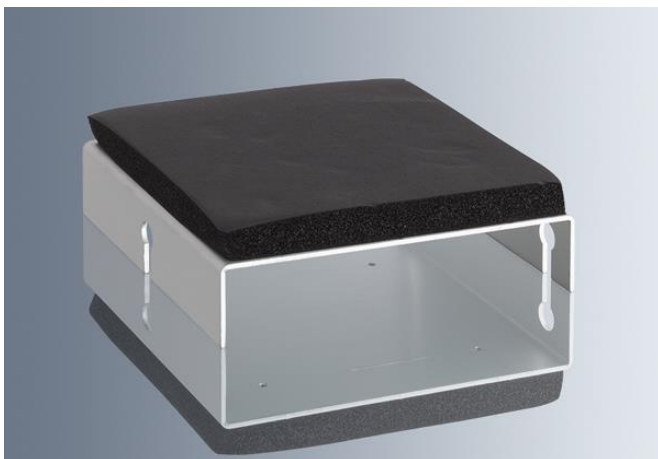**Single-hand shaking attachment
(Cat. No. 7204015):**

This attachment is suitable for shaking several test tubes or small round-bottom flasks up to 100 ml, while held in single hand - ideal for the simple and quick laboratory shaking need.

Recommended max. shaking speed:
1100 rpm

2. General information

The instrument is designed in accordance with safety class I. It is built and tested in accordance with DIN EN 61010. It left our warehouse working perfectly well and safely. To maintain the proper safety and operational function of the instrument, the user must respect the instructions and safety guidelines of this manual.

The instrument comes with a SCHUKO plug (DIN 49441, 16 A, 250 V). When operating the instrument in countries with different AC plug systems, use an approved adapter or have a qualified electrician replace the AC plug with an approved model suitable for the country of operation.

2.1 Unpacking the instrument

Unpack the instrument carefully and check to see that it is not damaged. It is important that any damage incurred during transport, such as indented sides of the carton, chemical impairment of the carton, water stains or other physical impacts which might have damaged the content, be recognized at the time of unpacking. Notify your carrier or forwarding agent immediately in case of such damage. You may unpack the carton carefully if there is no visible damage of the packing material.

2.2 Safety instructions

Note:

If the instrument is not damaged and all parts are complete, you may start to operate the device after reading the instruction manual.

Important note:

Read this instruction manual carefully and make sure everything is understood before operating the instrument. Should there be any further questions after reading the manual concerning the setup, operation or warranty, please contact:

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Attention:

Please comply with all safety and accident-prevention regulations applicable to laboratory work.

Caution:

When connecting the instrument to an AC power outlet, ensure that your local supply voltage corresponds to the specifications indicated on the instrument.

**Caution:**

The On/Off switch does not disconnect the instrument from the power source. Remove the plug from the AC power outlet to disconnect the instrument from the mains supply entirely.

**Attention:**

Do not open the instrument! Repairs are only to be carried out by trained service technicians. Any work on the electronics of the unit should only be done by knowledgeable and trained personnel. In the case of unauthorized operation as well as negligent or wilful damage, the warranty expires.

3. Intended use

The shaker is a multi-purpose shaker which can be used with several attachments for various kinds of shaking in chemical and pharmaceutical laboratories as well as in schools and universities.

To ensure maximum service life, observe the specified ambient conditions (temperature and humidity) and make sure that the instrument is not exposed to a corrosive atmosphere.

4. Setup and operation

Make sure to place the instrument on a level table. Allow for sufficient free space around the instrument to prevent anything from being knocked during the shaking operation.

Please observe the ambient conditions indicated in the technical data especially the voltage and observe the safety instructions.

**Caution:**

Make sure that there are no spills during shaking operation. Use test tubes which are high enough or reduce the fluid level or use a cover for the test tubes. For your own security wear protective clothing and glasses.

Place the attachment on top of the shaking unit and fasten the attachment by means of the three screws, one center screw on the front and two screws on the reverse side. All attachments are interchangeable and equipped with the same mounting system.

Switch on the instrument via toggle switch and choose the shaking function desired with the corresponding switch position:

- Switch position 0 (middle): The instrument is switched off.
- Switch position I (top): For continuous operation
- Switch position II (bottom): The instrument can be driven with time control. When choosing a time period by means of the left control button, both yellow lamps light up. The instrument starts shaking until time is up.

The required speed (continuously adjustable shaking frequency) can be set by means of the control button at the right side. Make sure that the shaking frequency is not set too high when starting to work. Otherwise there's a risk of collision with other objects or instruments placed next to the mixer. Always start with the lowest shaking frequency (100 rpm).

5. Cleaning and maintenance

The top of the shaking platform attachment may be cleaned with washing-up liquid and warm water. The adhesive mat can be easily cleaned with soap and water for long lasting use. It can be used again after drying.

The drive unit is driven by an eccentric shaking base by strong ball bearings and does not need service.

6. Dismantling and disposal

1. Switch off the instrument by means of the rocker switch.
2. Disconnect the instrument from the mains.
3. Now you can remove the instrument from the working area.

Damaged and/or disposed electric or electronic equipment has to be disposed of at the designated recycling depot. Also the packing material should be disposed of environment-friendly (material separation).

7. Transport and storage

7.1 Packing/Transport

Please use, if possible, the original packaging. Protect the instrument by means of bubble wrap against mechanical shocks from the outside. In case you do not use the original packaging, please mark the box with the following notes:

- Glass symbol (Fragile! Handle with care!)
- Umbrella (Keep dry!)
- Content (List of content)

7.2 Storage

For proper storage please observe the following conditions of surrounding:

max. ambient temperature: 5-40 °C
max. ambient humidity: 80 %

8. Warranty and Liability:

Warranty:

Paul Marienfeld GmbH & Co. KG provides a guarantee for a period of 24 months from date of delivery for the instrument purchased by you. This warranty applies only to defects in material or workmanship. There is no warranty for defects or malfunctions caused by neglect, improper use or improper service.

In case of defects in material or workmanship, the instrument will be repaired or replaced. In case of replacement, the warranty for the new unit will persist for the remaining period of primary warranty period. Charges for returning the instrument must be borne by the customer. Any claims regarding guarantee and liability are limited to professional and proper use of the instrument and observance of the general safety instructions and this instruction manual.

Liability:

Paul Marienfeld GmbH & Co. KG is not liable for any material or personal damages caused by improper use and does not accept liability for consequential damages. The warranty expires in case of removal or alteration of the identification label and/or the serial number of the instrument.

Attention:

In case of malfunction do not try to carry out any repair works. The instrument does not contain any part which may be serviced or maintained by the user. In case of any attempt by the user to repair the unit, the warranty expires!

Do not open the instrument. Any work on the electronics of the unit should be done by knowledgeable and trained personnel only!

9. Technical data

Power supply:	115 Volt or 230 Volt (see back panel)
Speed:	stepless adjustable from 100 to 1100 rpm
Timer:	from 2 to 60 minutes
Dimensions basic drive:	180 x 140 x 125 mm (L x W x H)
Weight basic drive:	approx. 4 kg
Input:	40 Watt
Admissible ambient temperature:	+5 °C up to +40 °C
Admissible max. humidity:	80 %
Maximum time limit:	100 %

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